# Levenhuk DTX TV LCD

## Digital Microscope

EN	User Manual
BG	Ръководство за потребителя
CZ	Návod k použití
DE	Bedienungsanleitung
ES	Guía del usuario
HU	Használati útmutató
IT	Guida all'utilizzo
PL	Instrukcja obsługi
PT	Manual do usuário
RU	Инструкция по эксплуатации

**220x**  
**ZOOOOM**

**3** MEGA  
PIXELS

**HDMI**  
**1080 P**

**levenhuk**  
Zoom&Joy


# Levenhuk DTX TV LCD

## Digital Microscope

EN

### Microscope parts

- |                                  |  |
|----------------------------------|--|
| 1 Up                             | 11 Focus wheel                             |
| 2 Down                           | 12 HDMI out                                |
| 3 Setup/Delete                   | 13 1/4" screw                              |
| 4 OK (Confirm)                   | 14 TV out                                  |
| 5 Power on/off                   | 15 DC: charge battery<br>USB: output to PC |
| 6 LED Dimmer                     | 16 Photo/Video switch                      |
| 7 Capture remote control unit in | 17 Capture                                 |
| 8 Charging indicator             | 18 Playback                                |
| 9 MicroSD slot                   |  |
| 10 2.4" color LCD display        |  |

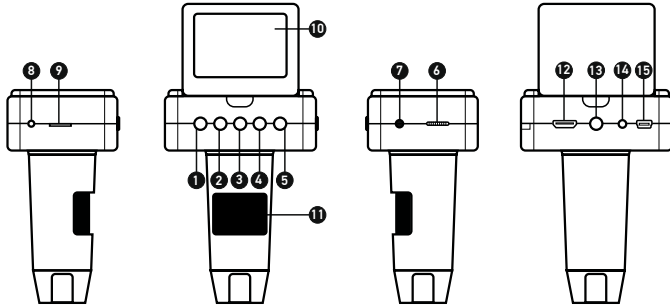
The kit includes: microscope, microscope stand, AC adapter, USB cable, HDMI cable, TV cable, capture remote control, 4G microSD memory card, software CD, user manual and warranty slip.

Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause **PERMANENT RETINAL DAMAGE** and may lead to **BLINDNESS**.

### Microscope assembly

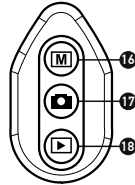Take the microscope and all of the accessories out of the package. Place the stage on a stable surface. Attach the mount locking ring (16) to the metal tube of the mount. Insert the metal tube into the opening on the stage and lock it in place with the ring (16). Insert the microscope into the clamp on the mount arm and tighten the locking screw (14). Tighten the locking knob (15) to lock the mount arm in place. To power the microscope connect it to your PC with a USB cable.

### Display icons

-  Capture an image
-  Start and stop video recording
- $\pm 0$  Exposure

08594 Total number of images that may be captured (based on the current resolution)

-  MicroSD status icon
- 3M Current image resolution
-  Battery status indicator

*Note: Interface icons will be hidden if no key has been pressed and held for 5 seconds.*

### PortableCaptureHD software

Insert the installation CD into your CD drive. The Main Menu window will be displayed. Click the **Install PortableCaptureHD** to install the application and follow the instructions on screen. The installation CD contains a copy of this user manual, which you can read by clicking the corresponding entry in the **Main Menu**. Connect the microscope to your PC with a USB cable and run the PortableCaptureHD application. A pop-up message will warn you if the microscope is not connected to the PC.

Capture an image

Start and stop video recording

Sequence capture (you can set the time to begin capturing images or videos, interval between them, video duration and the number of images or videos to be captured)

Full screen mode. To exit full screen mode, press **Escape** on your keyboard or double-click anywhere on the screen

## Using the microscope

### Connection to TV/monitor with HDMI input

- 1) Connect the microscope to the TV/monitor by inserting the HDMI cable into the HDMI input port. Turn on the power of your TV/monitor and select the HDMI setting.
- 2) Power the microscope with the supplied power adapter. If the microscope is not already on, turn on the power by pressing the Power on/off button (5), and then you can start the preview on the screen.

### Capturing an image

- 1) Insert a microSD card into the microSD slot (9). Maximum card size is 32GB.
- 2) Turn the Focus wheel (11) to focus the image of your object to its best focus.
- 3) Press the Capture button (17) on the capture remote control and the photo will be saved to the microSD card.

### Playback

- 1) Insert a microSD card in the slot.
- 2) Press the Playback button (18) on the capture remote control to switch to playback mode.
- 3) Press the Down button (2) to watch the previous captured photo. Press the Playback button (18) on the capture remote control again to return to preview.

### Deleting files

- 1) Insert a microSD card in the slot.
- 2) Press the Playback button (18) on the capture remote control to switch to playback mode.
- 3) Press the Setup button (3) to choose a file to delete.

### Mass storage on PC

- 1) Insert a microSD card in the slot.
- 2) Connect the microscope to your PC with the USB cable to download images from the card to your computer.

### Preview on PC

- 1) Only works with the software installed.
- 2) Remove the microSD card from the slot.
- 3) Connect the microscope to your PC with the USB cable.

## Display menu

Press Setup button (3) to open the main display menu. From this menu you can set image and video resolution, adjust image exposure, set the interface language, turn on the power saving mode, hide or display date and time, set up the TV connection, format the SD card, reset settings and check the hardware version. Also, this menu provides access to the Loop recording mode and Date Stamp option.

### Resolution

Select the image or video **Resolution** from the main menu and choose the required resolution by pressing the Down button (2), then confirm by pressing the OK button (4).

*Note: This model supports the following image resolutions: 3M, 5M, 8M, 12M; video resolutions: 1080P, 720P, VGA.*

### Exposure (Brightness)

Select **Exposure** from the main menu and then select a target value.

### Date Stamp

Select **Date Stamp** from the main menu and then select a target option.

### Date and time

Select **Date/Time** from the main menu by the Down button (2) and set the current date and time. Press the OK button (4) to confirm entry. You can also choose one of the following date formats: YY/MM/DD, DD/MM/YY or MM/DD/YY. Press the Setup button (3) to return to preview mode.

### Loop recording

Select the duration of the video clip to turn on the **Loop recording** mode. While the **Loop recording** mode is on, each video clip of selected duration will be recorded over the previous one.

### Power saving

Select **Auto Power Off** from the main menu and set the idle time before the microscope automatically shuts down. Press the OK button (4) to confirm selection.

### Language

Select **Language** from the main menu and choose one of the existing settings. Press the OK button (4) to confirm selection.

### TV connection

Select **TV Mode** from the main menu and then select, between NTSC and PAL, the correct mode compatible with your TV. Press the OK button (4) to confirm selection.

### MicroSD card formatting

Select **Format SD** from the main menu to format the installed memory card. **All the files on the card will be deleted after card formatting.** Normally card formatting is not necessary, unless your card is not able to store the captured image.

### Settings reset

Select **Default Setting** from the main menu to revert back to factory settings.

## Hardware version

Select **Version** from the main menu to check the current hardware version.

## Application menu

### File

**Photos Directory:** Set a directory for captured images.

**Videos Directory:** Set a directory for recorded videos.

### Options

**Resolution:** Set image resolution.

**Date/Time:** Hide or display date and time of capture during preview.

**Language:** Change the language of the user interface.

**XY Guide:** Hide or display a grid on images.

**Full screen mode:** Go full screen.

### Capture

**Photo:** Capture an image.

**Video:** Record a video.

### Capturing an image

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu.

### Recording a video

Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu.

### Sequence capture

Click the corresponding icon on the toolbar. A dialog window will appear where you can set a number of parameters for sequence capture.

Click the **Photo** radio button to set the time to begin capturing images, interval between them and the total number of images. Click the **Video** radio button to set the time to begin recording videos, their duration, interval between them and the total number of clips.

Save as

Copy to clipboard

Copy file

Previous image

Next image

Undo the last action

Repeat the last action

Draw

Text

Measurement

Calibration

## Calibration

Before calibrating the microscope, turn the grid display on. To do so, select **On** from **Options > XY Guide**. By default, the grid is visible. Place the calibration scale on the stage and focus your view. Make sure that the vertical axis of the grid is parallel to scale divisions and capture an image. Double-click the image thumbnail to open it in a separate window.

Click the corresponding icon on the toolbar to begin calibration. Using your mouse, choose two points on the image (you have to know the actual distance between these two points). Once you choose the second point, a dialog window will open, where you have to enter the known distance in the **Actual dimension** field. The application will automatically calculate the magnification of the image. Click **OK** to close the dialog window. The calculated magnification will be displayed in the **Magnification** field.

**Note:** Chosen points should form a horizontal line.

You can check calibration results by measuring the same distance with the **Straight line** tool. Click the measurements icon, choose the **Straight line** tool and draw a similar line on the image. If the measured distance is equal to the actual distance, calibration has been successful.

## Recalibration

The calibration process has to be repeated if the magnification or focus has been changed during observations.

Click the corresponding icon on the toolbar again and select **Reset picture magnification** from **Reset magnification**. Repeat the calibration process, as described above. Change the magnification and capture a few images. Double-click one of the thumbnails to open the resulting image in a preview window. Click the calibration icon on the toolbar and select **Set picture magnification** from **Set magnification**. Enter the current magnification in the dialog window that appears and click **OK**.

If you want to use the same magnification when capturing images in the future, click the calibration icon on the toolbar and select **Set capture magnification** from **Set magnification**.

## Measurements

It is recommended to calibrate the system before proceeding with measurements.

**Straight line.** Press and hold the left mouse button to start drawing a line. Release the left mouse button to complete the line.

**Continuous Line.** Press and hold the left mouse button and draw a line. Release the left mouse button to complete the segment. You can continue adding more segments to the free-form line.

**Radius Circle.** Draw a straight line on the image. The application will automatically calculate the radius, as well as circumference and area of the corresponding circle.

**Diameter Circle.** Draw a straight line on the image. The application will automatically calculate the diameter, as well as circumference and area of the corresponding circle.

**Three Points Angle.** Press and hold the left mouse button to draw a line on the image. Release the left mouse button and draw another line. The application will automatically calculate value of the resulting angle.

## Draw and text

PortableCaptureHD allows you to add drawings and textboxes to your images. Open an image in a preview window and click the **Pencil** icon on the toolbar. Select one of the tools from the drop-down menu and draw something on the image. By clicking the **TT** icon on the toolbar, you can edit font type and color to be used in textboxes.

## Working with the application for Mac OS users

Browse to the **mac** folder on the installation CD and copy **PortableCaptureHD** file to the desktop. Use a USB cable to connect the microscope to your computer. Run the PortableCaptureHD application by double-clicking its icon.

A magnified image of the observed specimen may be seen in the main window of the application. From the **Options** menu, select **Preview Size** to set the image resolution. Keep in mind that the preview window cannot be resized (standard resolution is 640x480). The image resolution is shown in the bottom left corner of the preview window.

From the **Options** menu, select **Date/Time** to display or hide the date and time of capture during preview.

Click the corresponding icon on the toolbar or select **Photo** from the **Capture** menu to capture an image. An image thumbnail will appear on the right side of the main window. Click the corresponding icon on the toolbar or select **Video** from the **Capture** menu to start recording a video. A video thumbnail will appear on the right side of the main window.

***Note:** During recording, the camcorder icon on the toolbar will turn red. Click the icon again to stop recording.*

Image and video thumbnails are displayed on the right side of the main window of the application.

Double-click an image thumbnail to open it in a separate preview window.

Editing images with PortableCaptureHD application is similar to working with other applications in Mac OS.

From the main menu, select **Quit PortableCaptureHD** to exit the application.

## Care and maintenance

**Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.** White LEDs are very bright; do not look at them as this may cause permanent damage to your vision. Take necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions. After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection. Do not try to disassemble the device on your own for any reason. For repairs and cleaning of any kind, please contact your local specialized service center. Protect the device from sudden impact and excessive mechanical force. Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws. Do not touch the optical surfaces with your fingers. To clean the device exterior, use only special cleaning wipes and special optics cleaning tools from Levenhuk. Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush. Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity. Be careful during your observations, always replace the dust cover after you are finished with observations to protect the device from dust and stains. If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope. Store the device in a dry, cool place away from hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures. To avoid data loss or equipment damage, always close the application first, and then disconnect the microscope from your PC. Wait until the illumination goes off completely, and disconnect the microscope. **Seek medical advice immediately if a small part or a battery is swallowed.**

## Battery safety instructions

Always purchase the correct size and grade of battery most suitable for the intended use. Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types. Clean the battery contacts and also those of the device prior to battery installation. Make sure the batteries are installed correctly with regard to polarity (+ and -). Remove batteries from equipment that is not to be used for an extended period of time. Remove used batteries promptly. Never attempt to recharge primary batteries as this may cause leakage, fire, or explosion. Never short-circuit batteries as this may lead to high temperatures, leakage, or explosion. Never heat batteries in order to revive them. Remember to switch off devices after use. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning.

## Specifications

Optics material	glass
Megapixels	3
Magnification	10–220x
Available image resolutions	3M, 5M, 8M, 12M
Available video resolutions	1080P Full HD, 720P, VGA
Photo/video format	*.jpg/*.avi
Focusing	manual, 10–50mm
Illumination	8-LED system with variable brightness
Power source	built-in Li-ion battery or 5V DC-in
Software language	English, German, Spanish, Russian, Italian, French, Portuguese, Dutch, Polish, Japanese, Korean, Chinese
Ability to connect additional equipment	support of microSD cards with capacity up to 32GB
Operating temperature range	–5...+50 °C / 32...122 °F

Levenhuk reserves the right to modify or discontinue any product without prior notice.

## System requirements

Operating system: Windows 7/8/10, Mac 10.7 and above

CPU: at least P4 1.8GHz or above, RAM: 512MB, video card: 64MB

Interface: USB 2.0, CD drive, any monitor with HDMI input.

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime warranty** against defects in materials and workmanship.

Lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from date of retail purchase. The warranty entitles you to free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all warranty conditions are met. For further details please visit our web site:

[www.levenhuk.com/warranty](http://www.levenhuk.com/warranty)

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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pericolosi e altri prodotti chimici, da apparecchi di riscaldamento, da fiamme libere e da altre fonti di calore. When using the microscope, try not to use it near flammable materials or substances (benzene, paper, cardboard, plastic, etc.), as the base may heat up during use, and might become a fire hazard. Always unplug the microscope from a power source before opening the base or changing the illumination lamp. Regardless of the lamp type (halogen or incandescent), give it some time to cool down before trying to change it, and always change it to a lamp of the same type. Always use the power supply with the proper voltage, i.e. indicated in the specifications of your new microscope. Plugging the instrument into a different power outlet may damage the electric circuitry of the microscope, burn out the lamp, or even cause a short circuit. **In caso di ingestione di una parte di piccole dimensioni o di una batteria, richiedere immediatamente assistenza medica.**

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lamp. Regardless of the lamp type (halogen or incandescent), give it some time to cool down before trying to change it, and always change it to a lamp of the same type. Always use the power supply with the proper voltage, i.e. indicated in the specifications of your new microscope. Plugging the instrument into a different power outlet may damage the electric circuitry of the microscope, burn out the lamp, or even cause a short circuit. **Procure um médico imediatamente se uma peça pequena ou uma pilha for engolida.**

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Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006—2021 Levenhuk, Inc. All rights reserved.